

ACTIGRAPH/ACTILIFE QUICK START GUIDE

This guide has been created by Arizona Performance Lab (apl.arizona.edu) under the guidance of Dr. Kristen Renner, and was adopted and edited by the Sensor Lab (sensorlab.arizona.edu) under the guidance of Dr. Gustavo Almeida. If you have any questions or edits to the guide please email Gustavo Almeida at gustavo@arizona.edu

Installation on a New Device

1. Go to this website to download the latest version of ActiLife 6
<https://actigraphcorp.com/support/software/actilife/>
 - . This website includes info on correct installation steps
<https://actigraphcorp.force.com/support/s/article/Installing-Reinstalling-ActiLife-5-x-6-x>
 - b. Product Key for Full Software: [Contact us](#)
 - c. Product Key for Lite Software: [Contact us](#)
 - . This key can be used in up to 5 devices
 - ii. If we want to move installation of ActiLife from one computer to another go here
<https://actigraphcorp.force.com/support/s/article/Installing-Reinstalling-ActiLife-5-x-6-x>
 - d. Online Tutorials <http://actigraphcorp.com/support/tutorials/>
 - e. General Manual <http://support.theactigraph.com/dl/ActiLife-manual>
 - f. Manual for ActiGraph GT9X Link + ActiLife
https://s3.amazonaws.com/actigraphcorp.com/wp-content/uploads/2020/03/05155628/ActiGraph_Link_UserGuide_E.200.6001_Revision6_FINAL.pdf
 - g. After you have correctly entered your Product Key and clicked on Activate, you will be prompted to enter your Name, Email, Organization, and Phone Number to complete the registration process when you open it for the first time. All of these fields are required.
2. You should eventually see this page


ActiLife v6.13.4 Lite Edition - No Devices Connected

File Edit Tools Help

Devices Wear Time Validation Scoring Sleep Batch Sleep PLM Graphing NHANES GPS Feature Extraction CentrePoint Data Vault

Initialize Download Refresh Refresh All Identify Advanced

Automatic Refresh

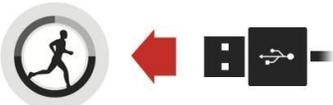


 **Introducing CentrePoint Study Admin**
Cloud-Based Study Management & Real-time Data Monitoring

The CentrePoint Study Admin system is ActiGraph's secure cloud-based technology platform designed to streamline workflows, simplify operational tasks, and solve logistical challenges associated with large, multi-site research studies and clinical trials.

[Learn more](#)

Initialization

1. Open the ActiLife software
2. Connect the ActiGraph dock to the PC using USB cable
3. Insert Link into dock with logo facing up – once it is connected it should appear in the grid on ActiLife under the ‘Devices’ tab
4. Select ‘Initialize’ and then select ‘Regular Initialization’
 - a. The most used initialization parameters are pre-selected as system defaults
 - b. You should have 30 Hz as the sampling rate, Enable IMU selected, and the start time set to be around the time your subject will arrive and have the stop time be set for around 2-3 hours after depending on the length of the study to be safe.
 - c. Usually after the GT9X Link is reinitialized the time display would be disabled and the screen is blank. Make sure to check ‘**Show Display**’ before moving forward.
 - d. Enter the subject info with their subject name, gender, height and weight if you know it or have the subject there, their date of birth, and their non-dominant wrist
 - i. The subject name is required for initialization, but the other information can be entered during download or data analysis

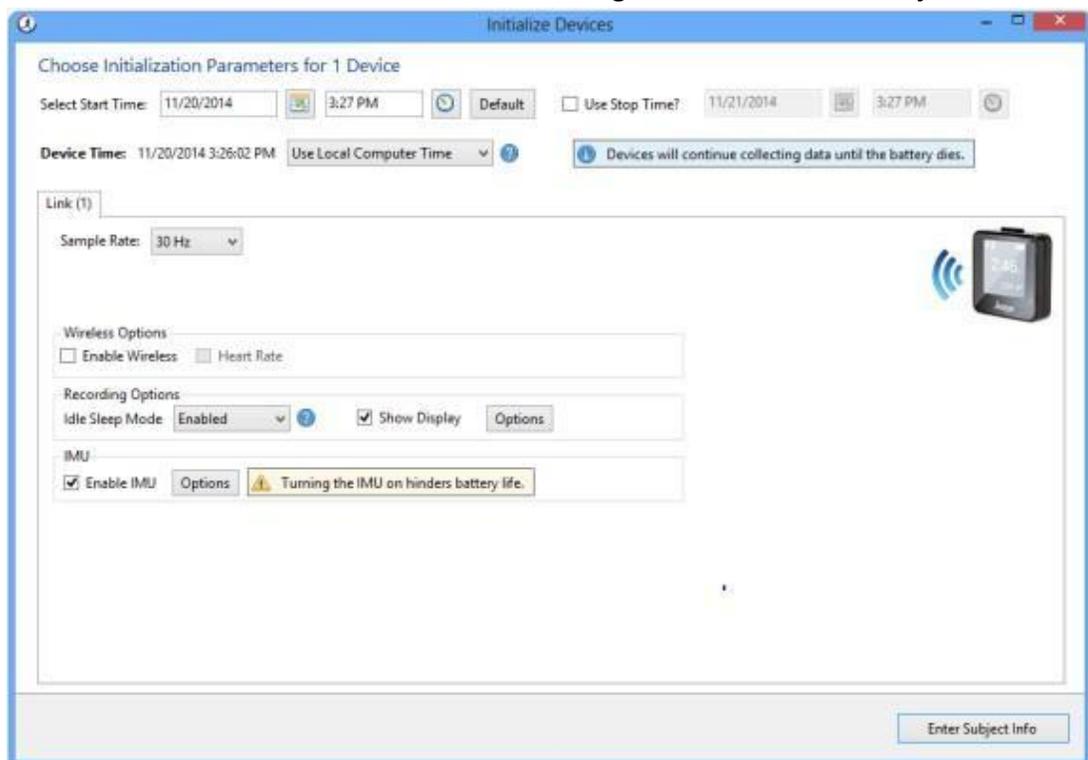


5. Select ‘Initialize 1 Device’. A progress bar in the device grid will indicated when the initialization process is completed
6. After initialization is complete, remove the ActiGraph Link from the dock. Once removed, a 10 second wear sensor calibration countdown will begin. For the watch

immediately insert the device into the watch band and lay it down on a flat surface until the countdown is complete

Downloading Data

1. Open the ActiLife software
2. Connect the ActiGraph dock to the PC using USB cable
3. Insert Link into dock with logo facing up – once it is connected it should appear in the grid on ActiLife under the ‘Devices’ tab
4. Verify the box in front of the device is checked and select ‘Download’ from the taskbar menu
5. A dialog box will open to display the download options listed below. The most commonly used download parameters are preselected as defaults

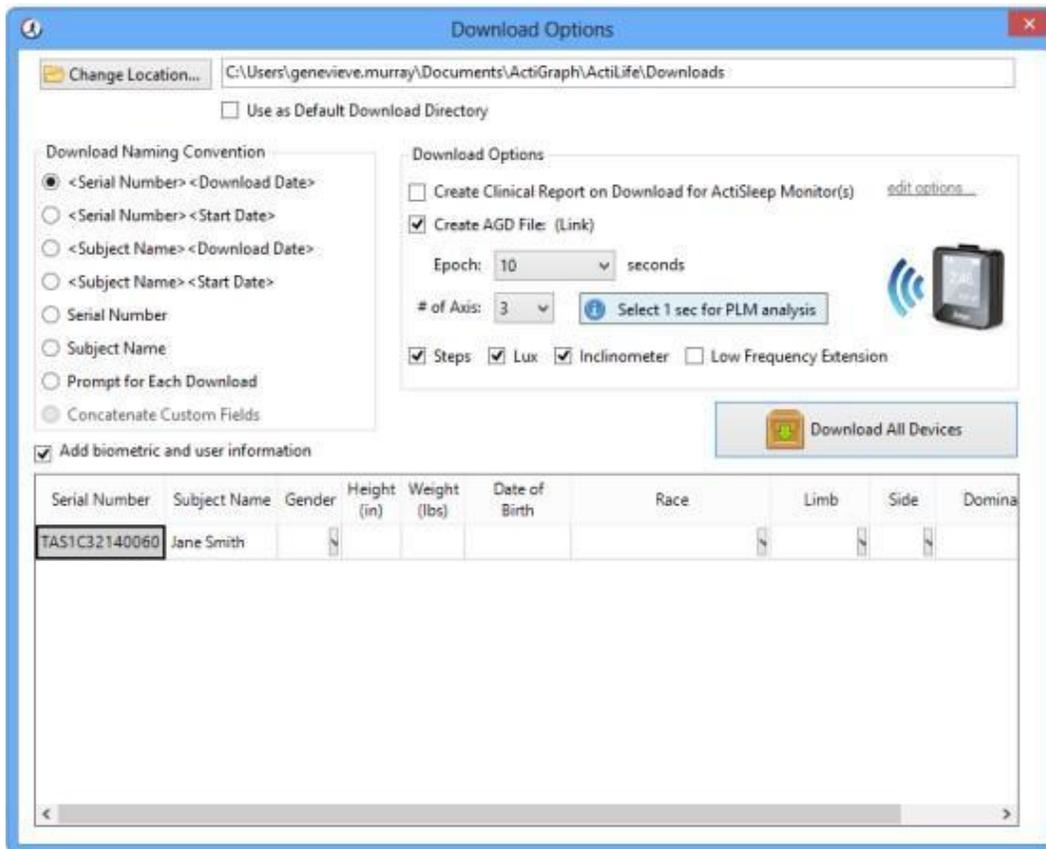

Serial Number	Subject Name	Gender	Height (in)	Weight (lbs)	Date of Birth	Race	Limb	Side	Dominance
TAS1C32140060	Jane Smith	[dropdown]	[dropdown]	[dropdown]	[dropdown]	[dropdown]	[dropdown]	[dropdown]	[dropdown]

6. Once Download Options are selected, click ‘Download All Devices’. The dialog box will close and a progress bar will appear under the ‘Status’ column in the grid
7. When the download is complete, a ‘finished downloading’ link will appear. Click this link to reveal options to view data, export the raw data file, and navigate to the download folder
8. After completing the download, select File->Import/Export/Convert->IMU in the ActiLife taskbar
9. Select the raw (.gt3x) file from the list and click ‘Open’

10. A status bar will appear while ActiLife converts the raw file to CSV. Once completed, the file location window will open and the IMU file will appear in the list
11. Once you have everything downloaded you will upload it to the Box and should have a 60sec.csv, 60sec.dat, IMU.csv, .gt3x, 10sec.agd, 60sec.agd, 10sec.csv, and 10sec.dat file

ActiLife Software

Support Center / Software / ActiLife Software



Software

Current Release: v6.13.4
Release Date: Apr 22, 2019

Download ActiLife 6

Related Downloads

- 9 [Net 4.0 Full \(Online Installer\)](#)
- 9 [Net 4.0 Full \(Offline Installer\)](#)

Release Notes

Version 6.13.4 – April 2019

CT3X file format support for the CentrePoint Insign Watch (CPW)

Version 6.13.3 – August 2016


This release addresses several usability issues with initialization and the API.

- Show busy dialog when a device disappears during an automatic firmware update.
- Store wear location for all devices that can accept it, not just Link.
- Show product key when deactivating without an internet connection.
- If a user's DF setting is over 100%, ActiLife scales correctly.
- Link initialization with subject feedback should always enable epoch collection.
- Ensure the API doesn't hang during automatic firmware updates.
- Calling USBList in the API works when there are more than one device connected.
- ActiLife no longer crashes when creating a clinical report if there are no days with wear time.
- In a shared user environment, ActiLife can now detect if it is already running on another user's login without crashing.
- When loading sleep periods for an ACD, ignore ones that aren't in the file's epoch range.

Retrieving Your Version of ActiLife

If your product key requires you to run a certain version of ActiLife, you can find it here. Provide just your product key in the new page and you'll be given a link to the version of ActiLife that your product key is able to be used with.

[Retrieve ActiLife Installer by Product Key](#)

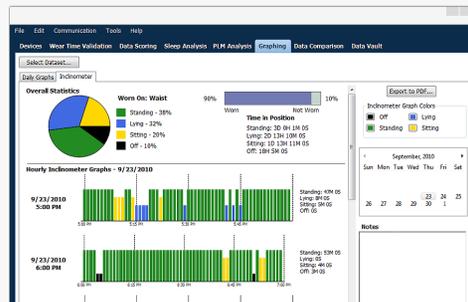



Solutions

- Clinical Trials
- Academic Research
- Sleep Assessment

Products

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
- CP Insight Watch
- wGT3X-BT
- CentrePoint
- ActiLife

Company

- About Us
- News & Events
- Leadership
- Contact Us

Support

- Support Center
- Downloads
- Tutorials
- RMAs



Copyright © 2018 ActiGraph Corporation. All rights reserved. [Terms of Use](#) | [Privacy Policy](#) | [Help](#) | [Contact Us](#)

